

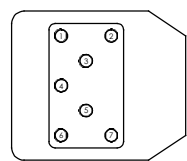


AMPERAGE DRAW PER LIGHT FIXTURE:
0.75A @ 230VAC

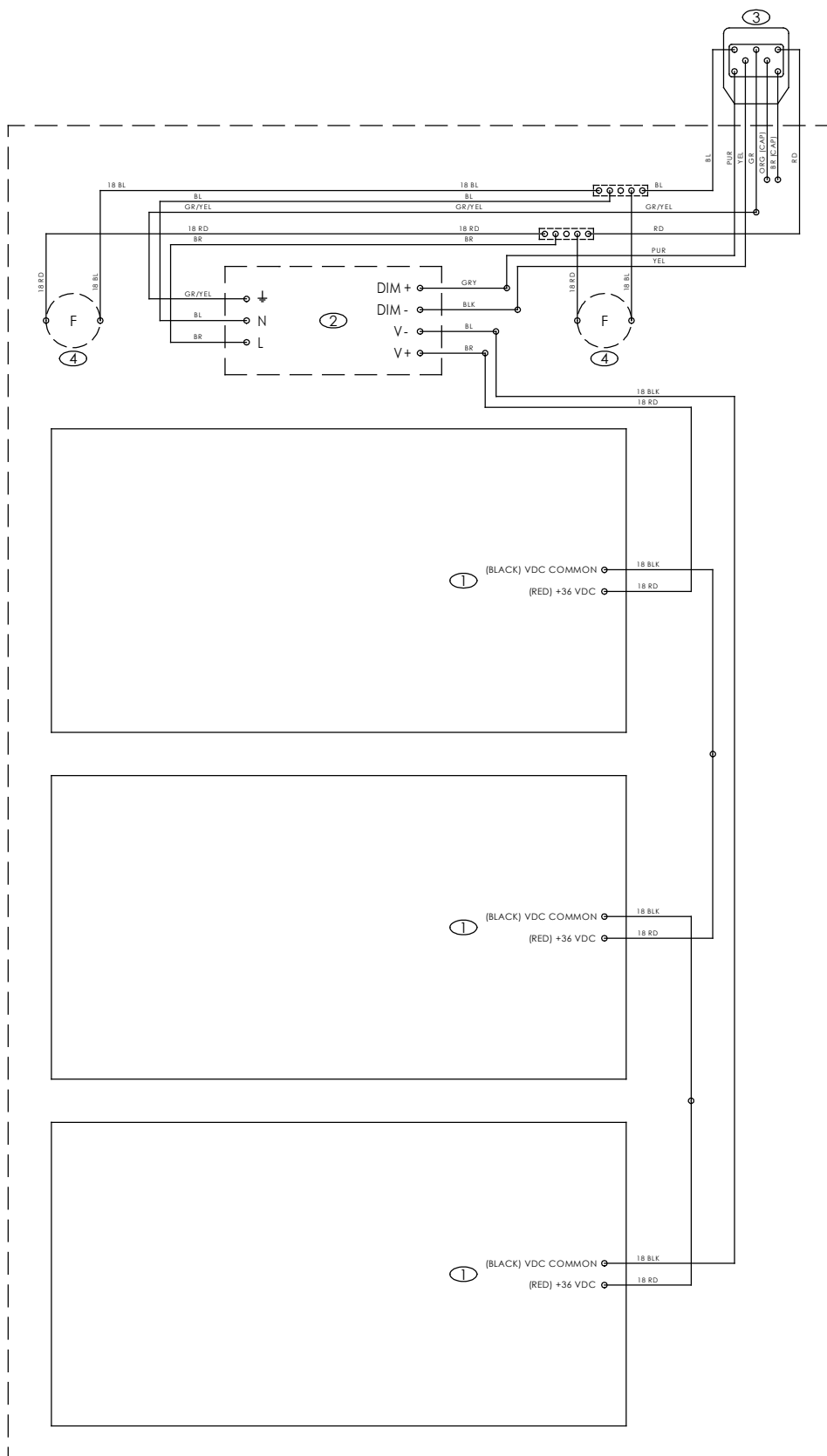

PIN 1 - RED
PIN 2 - BROWN
PIN 3 - ORANGE
PIN 4 - GREEN
PIN 5 - YELLOW
PIN 6 - BLUE
PIN 7 - PURPLE

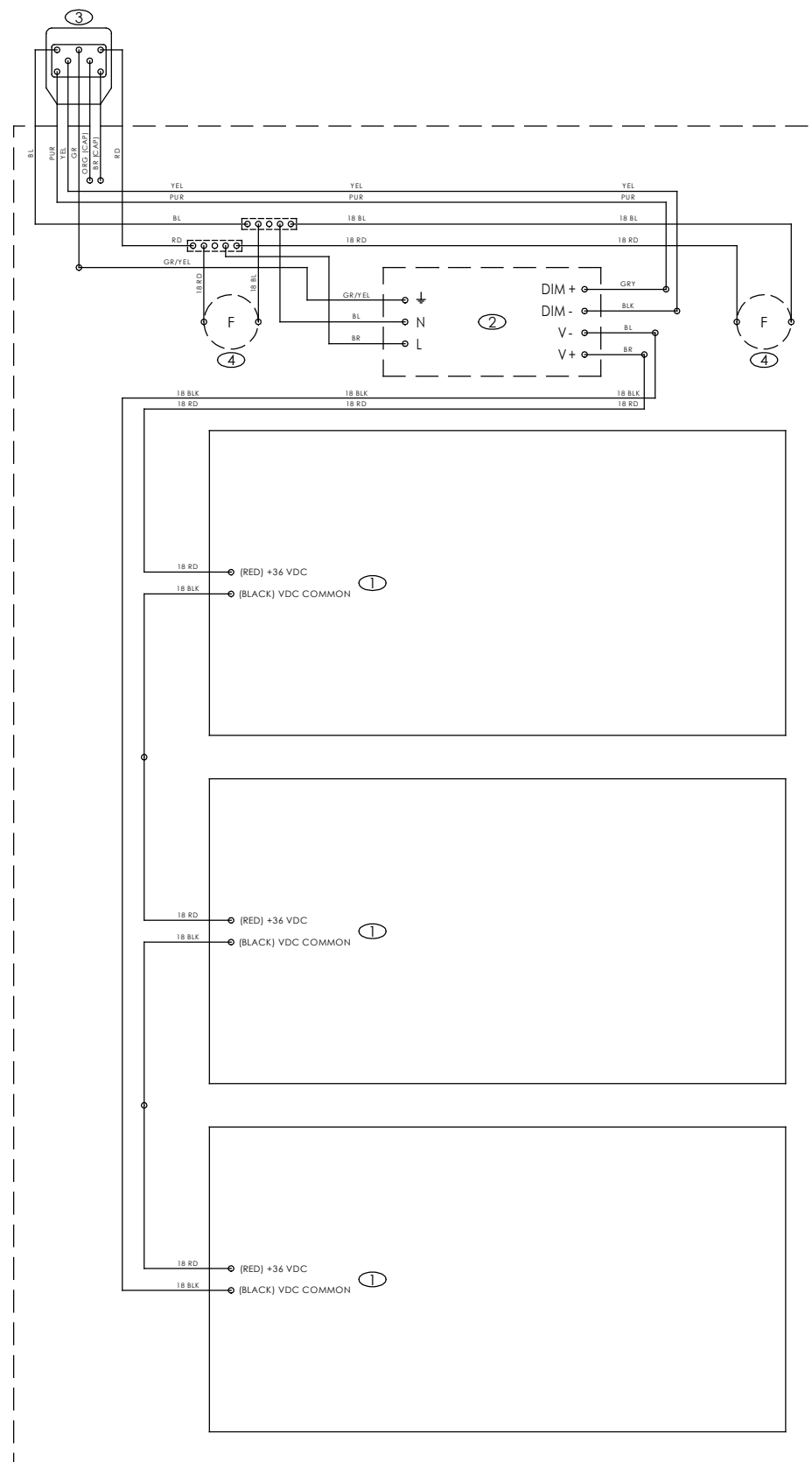

NOTES
1. CORD LENGTH TO BE 18".

- 1 - LED MODULE - DIMMABLE, WHITE, 36 VDC (ELD-066)
- 2 - LED DRIVER - HLG-185H-C1400B (ELD-071)
- 3 - PLUG - EMF #P-7-1, 7-PIN, 7-COLOR (EPL-052)
- 4 - LAMP BANK COOLING FANS - AXIAL, 53 CFM, 230VAC, 0.06A (EFA-111)

TITLE: ELECTRICAL SCHEMATIC - GENERAL LIGHTING			
MODEL: LT-36VL	DATE: 01/29/18	DRAWING: 2018-4-0-2	REV: B
BY: LCB			