


## IntellusUltra Key Definitions

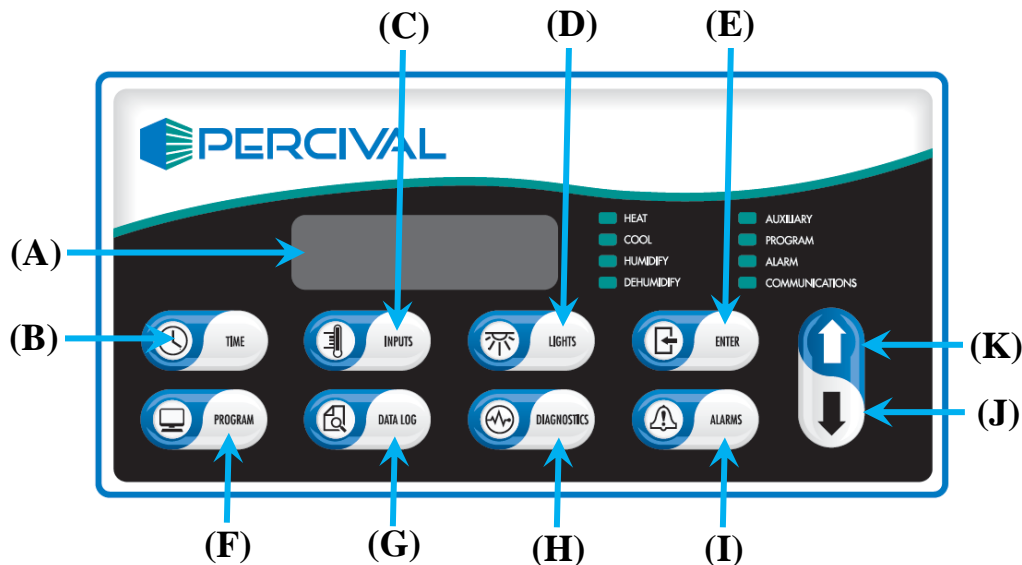

**Main Display (A)** - The main display shows all status and programming information.

**Time Key (B)** - The Time key is used to set the system time.

**Inputs Key (C)** - Used to set manual settings for temperature and humidity.

**Lights Key (D)** - This key is used to set lighting settings for the SciWhite® / PetriClear® LED tiles.

**Enter Key (E)** - This key is used to begin editing values and to accept changes of values.

**Program Key (F)** - This key is used to create/edit programs and to run programs or manual settings.

**Data Log Key (G)** - This key does not function on the standard IntellusUltra.

**Diagnostics Key (H)** - This key is used to enter the Diagnostics menu which allows you to view diagnostic information for the chamber.

**Alarms Key (I)** - This key is used to enter the Alarms menu. The Alarms menu contains entries for setting high and low alarm settings for temperature (Limit 1) and the redundant temperature limit (Limit 2).

**Decrement (J) / Increment (K) Keys** - These keys are used to scroll through available menu options and to decrease or increase values.